


Add or Remove Tests From Order

Manually Removing a Test from Order During LIS Downtime

1. Once logged into the **cobas® infinity** software, select the **Main** tab
2. Click on **Menu**
3. Navigate to **Order entry**
4. Select **Order Entry**
5. Type in existing **Order ID** to bring up patient
6. Double click on the test you wish to remove at bottom of screen

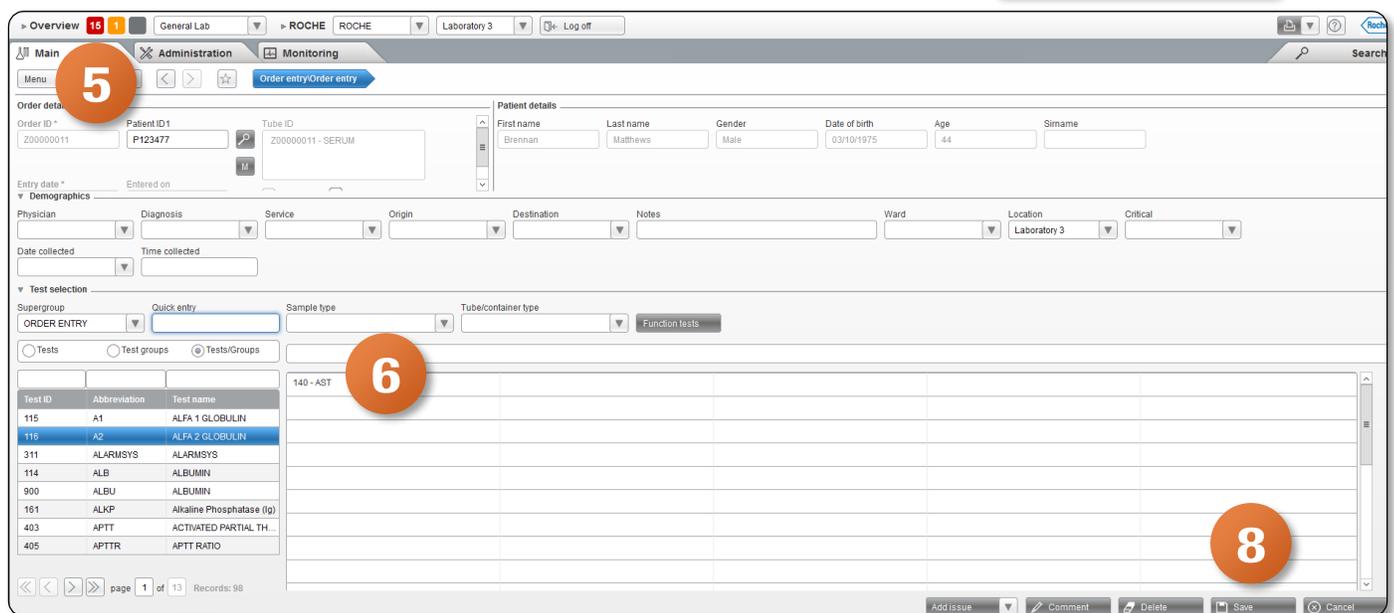


7. Click **Yes** to confirm you really want to delete the selected test
8. Click on **Save** to update the order



Add or Remove Tests From Order

Manually Adding a Test from Order During LIS Downtime

1. Once logged into the **cobas® infinity** software, select the **Main** tab
2. Click on **Menu**
3. Navigate to **Order entry**
4. Select **Order Entry**
5. Type in existing **Order ID** to bring up patient
6. Select a test to add to the order by double-clicking
7. Click on **Save**

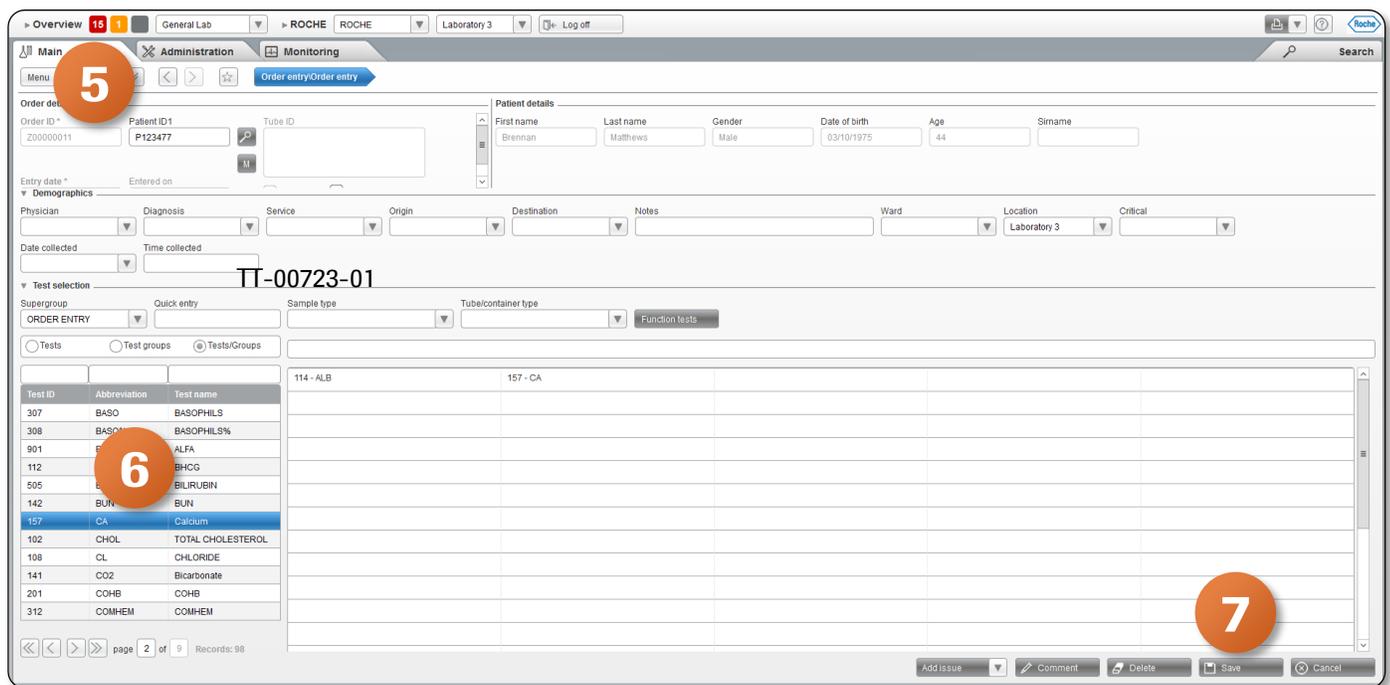

8. Once saved, **Order ID** and **Patient ID** will go blank indicating the system is ready for another entry

