


Connect, Login, and Logoff

Connect and Login to cobas® infinity laboratory solution


1. Open up desired Internet Browser. The **cobas® infinity** software is compatible with Chrome, Firefox and Internet Explorer
2. Type in the IP address of the server. Then add /infinity. We suggest making this a shortcut so it is saved for easy navigation in the future

Please Note: Your IP address will be different than what we are using in this example

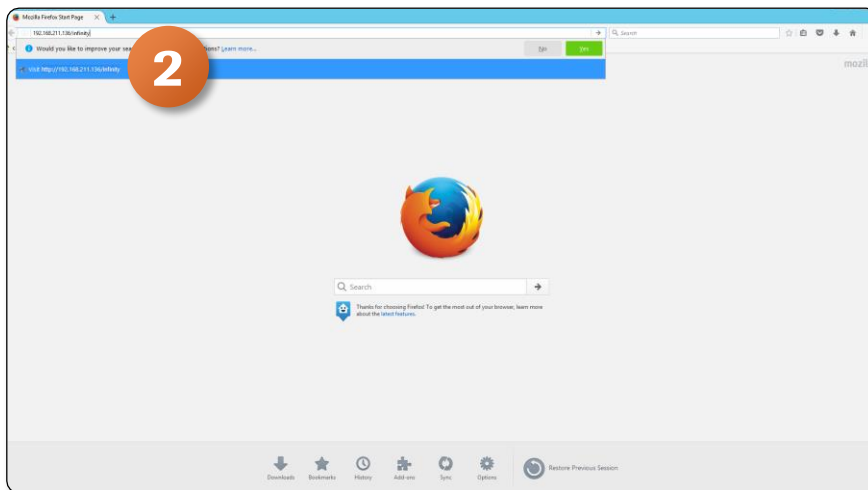

3. The **cobas® infinity** software Log on screen will now appear
4. Type in the **User name** and **Password**
5. Click on **Confirm**

Please note: If you exceed the maximum allowed attempts to login, you will need to contact your administrator to reset your password.

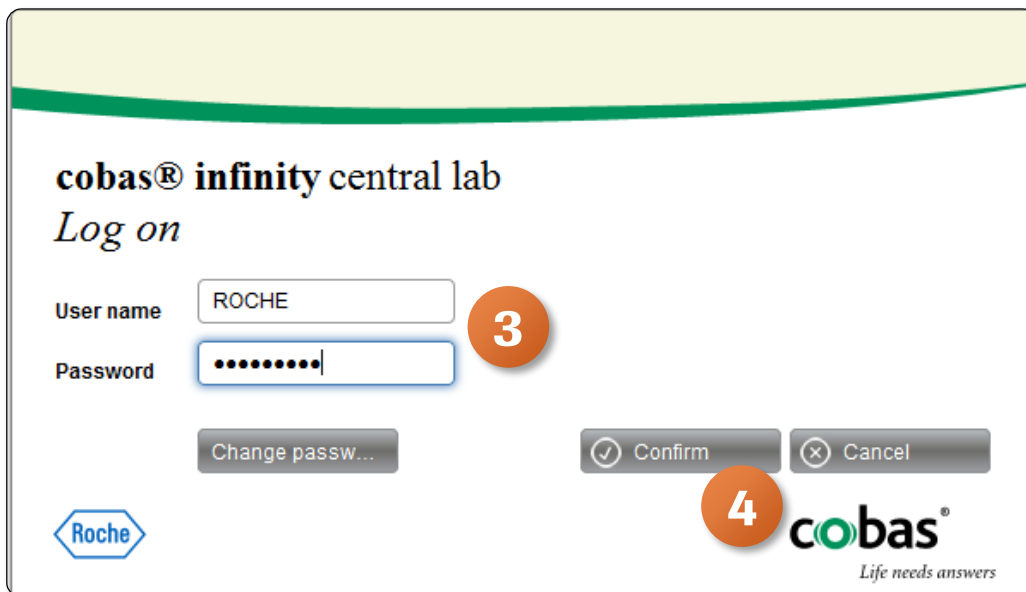

Connect, Login, and Logoff

Logoff of cobas® infinity laboratory solution

1. From any screen, choose the **Log off** button located in the Overview bar
2. To confirm that you want to exit the software, select **Yes** from the callout

