


Exporting Information

1. Once logged into the **cobas[®] infinity** software, navigate to desired screen
2. From the Overview menu bar, select the drop-down list
3. From the Grid group box, choose one of the following options:
 - Export to TXT: produce a text file using the current browser
 - Export to HTML: produce a web file using the current browser
 - Export to PDF in Vertical or Horizontal format: produce a PDF file using the current PDF viewer

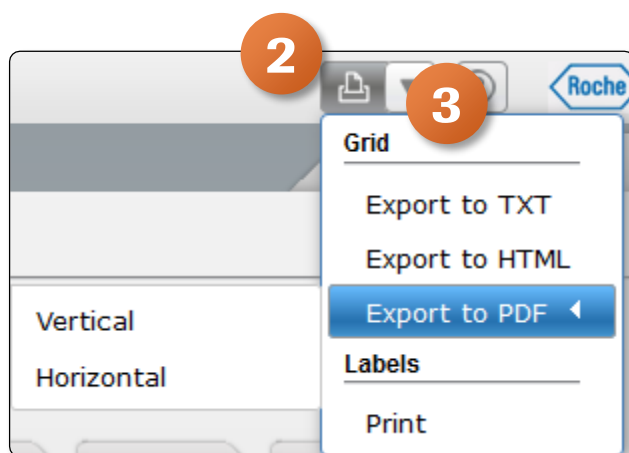

4. Once the report is generated, right click to see options. Select **Save Page As** to save this document
5. The button performs the last selected action. To print tables again, click print button or press F5

Screen report
Created by ROCHE
Date/Time of query 08/27/2019

Priority	Entry Date	SampleID	Entry Tim	Last name	First name	Patient ID	WA Status	Chemistry	Urines	# Tests	TAT	Critical	Alarms
General La	06/21/2019	Z00000011	19:41:42	Matthews	Brennan	P123477	2	1	1	7	-		0
General La	07/02/2019	Z00000012	09:30:41	Matthews	Brennan	P123477	2	1	1	1	-		0
General La	07/02/2019	Z00000014	09:39:01	Douglas	Estrella	P123480	2	2	1	11	-		0
General La	07/11/2019	Z00000015	09:31:54	Smith	Vern	000008	2	1	1	1	-		0
General La	07/11/2019	Z00000016	10:54:45	Smith	Adam	000009	2	2	1	11	-		0
General La	07/18/2019	Z00000021	10:20:37	Smith	Gene	000014	2	2	1	11	-		0
General La	07/18/2019	Z00000040	10:32:40	Smith	Adam	000009	2	2	1	2	-		0
General La	07/26/2019	Z00000030	13:24:55	Morrow	Anne	P123473	2	2	1	16	-		0

4 Right-click context menu options: Save Page As..., View Background Image, Select All, View Page Source, View Page Info, Inspect Element (Q)