
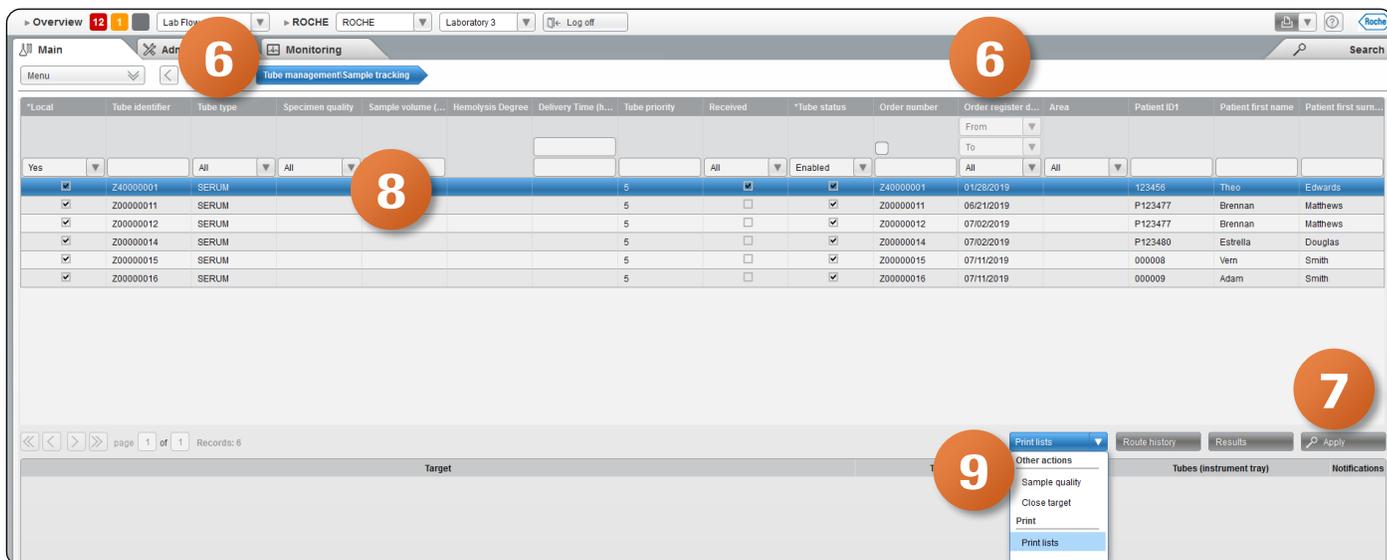


Indicating Sample Quality

If there is a problem and the sample cannot be used, it is useful to mark the quality of the sample, e.g. lipemic, hemolyzed, as correct.

1. Once logged into the **cobas[®] infinity** software, confirm that you are **Lab Flow** application
2. Select the **Main** Tab
3. Click on **Menu**
4. Navigate to **Tube management**
5. Select **Sample tracking**
6. Enter the Tube type and the Order registered
7. Click on **Apply**
8. Select the tube whose sample quality you want to indicate
9. Choose **Sample quality** from the drop-down menu

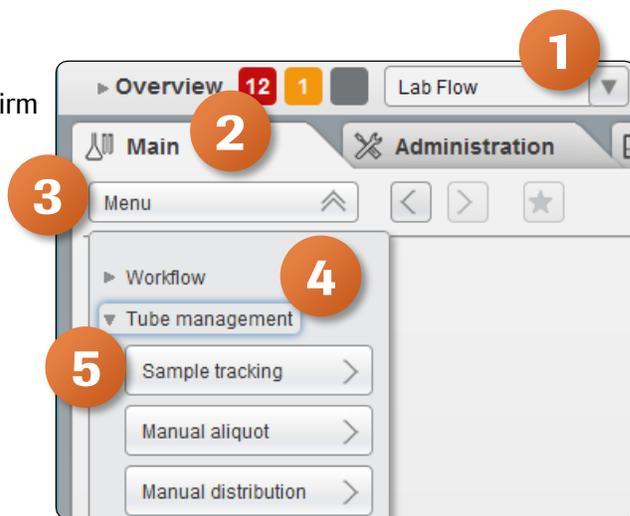

10. From the callout, select the **Sample quality** drop-down menu, select the desired option
11. Click on **Confirm**

