


Modifying Monitoring Screen

To begin you will need to figure out the exact name of the Work area you will be modifying

1. Once logged into the **cobas[®] infinity** software, select the **Main** tab
2. Click on **Menu**
3. Navigate to **Work areas**
4. This will show you a list of the Work area's for your laboratory

Please note: The Work area names are customized per laboratory What you see in our example won't necessarily reflect the names of your Work areas. The above navigation will help you to see the names of all Work areas you have created.

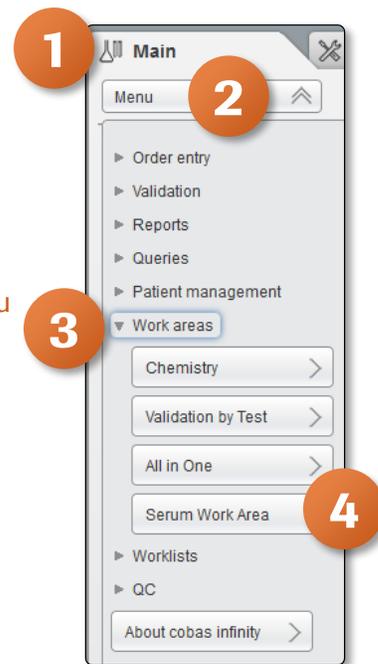


Once the Work area name has been confirmed, the next step is to confirm the exact name of the Monitoring screen

5. Navigate to the **Administration** tab
6. Click on **Menu**
7. Navigate to **Work areas**
8. Select **Entries**




Modifying Monitoring Screen

- 9. Double click on **Serum Work Area**
- 10. This populates the bottom portion of the screen. Click on the **Monitoring** tab

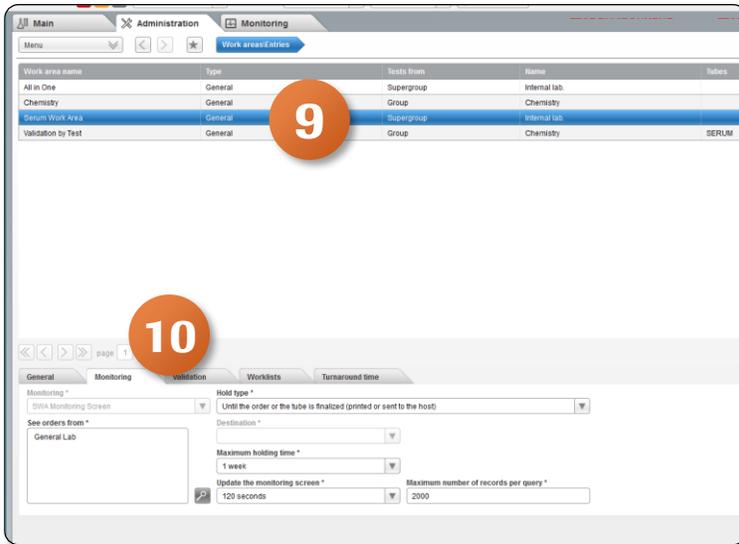

- 11. The first box in the Monitoring tab shows you what screen is associated with that Work area. In this example it is **SWA Monitoring Screen**

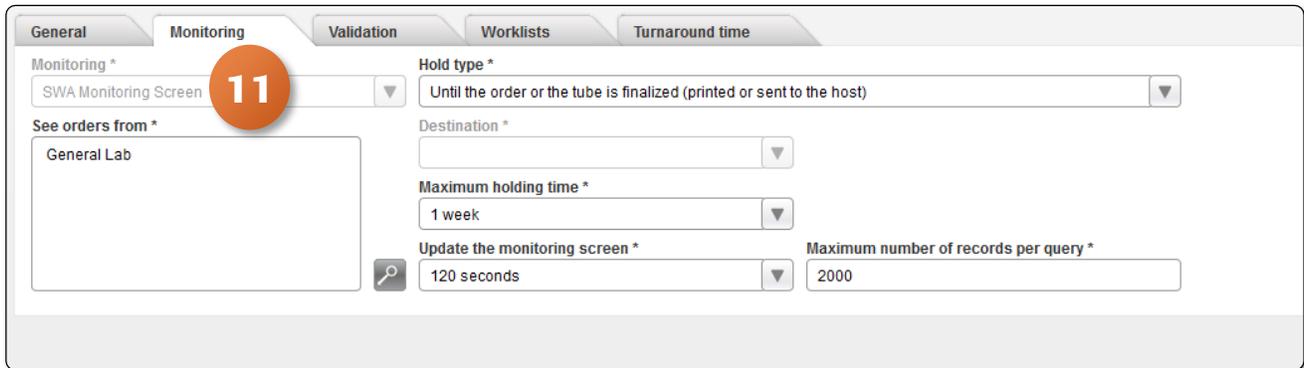


Modifying Monitoring Screen


Editing the Monitoring Screen

Now that we have the exact names of the screen, we can begin modification

1. Navigate to the **Administration** tab
2. Click on **Menu**
3. Select **General**
4. Click on **Configurable screens**

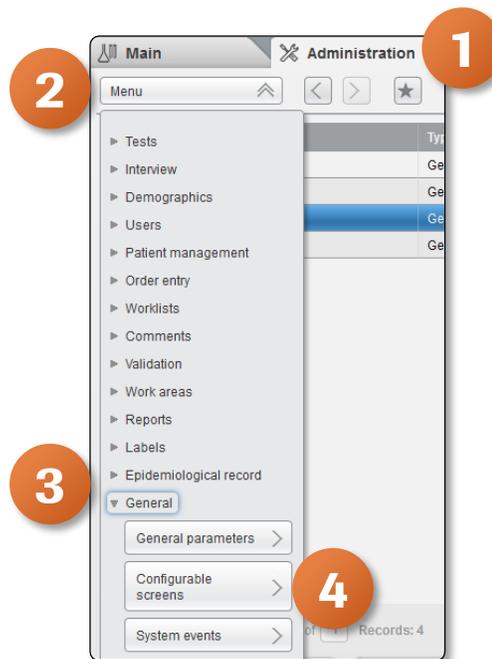


5. Select the Monitoring Screen associated with your Work area
6. You may need to use the arrows to see more options
7. Click on **Configure**

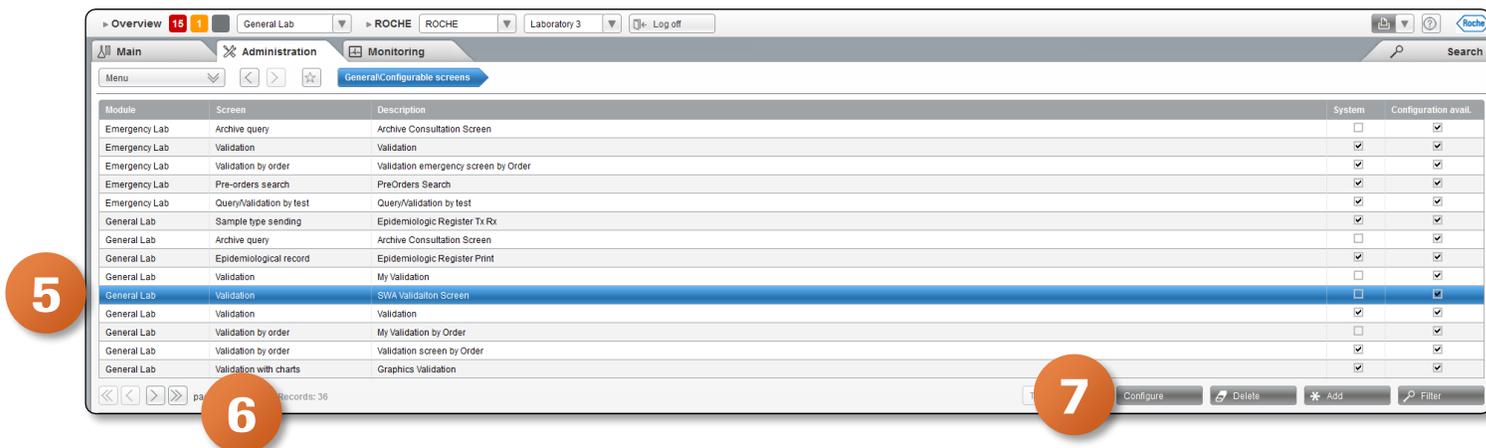

Modifying Monitoring Screen

- Click on **Add** to customize the Monitoring Screen
- Select the **Field type** and **Label**
- You have the option to change the **Width** of the columns by adding or reducing the number
- Click on **Confirm** once complete

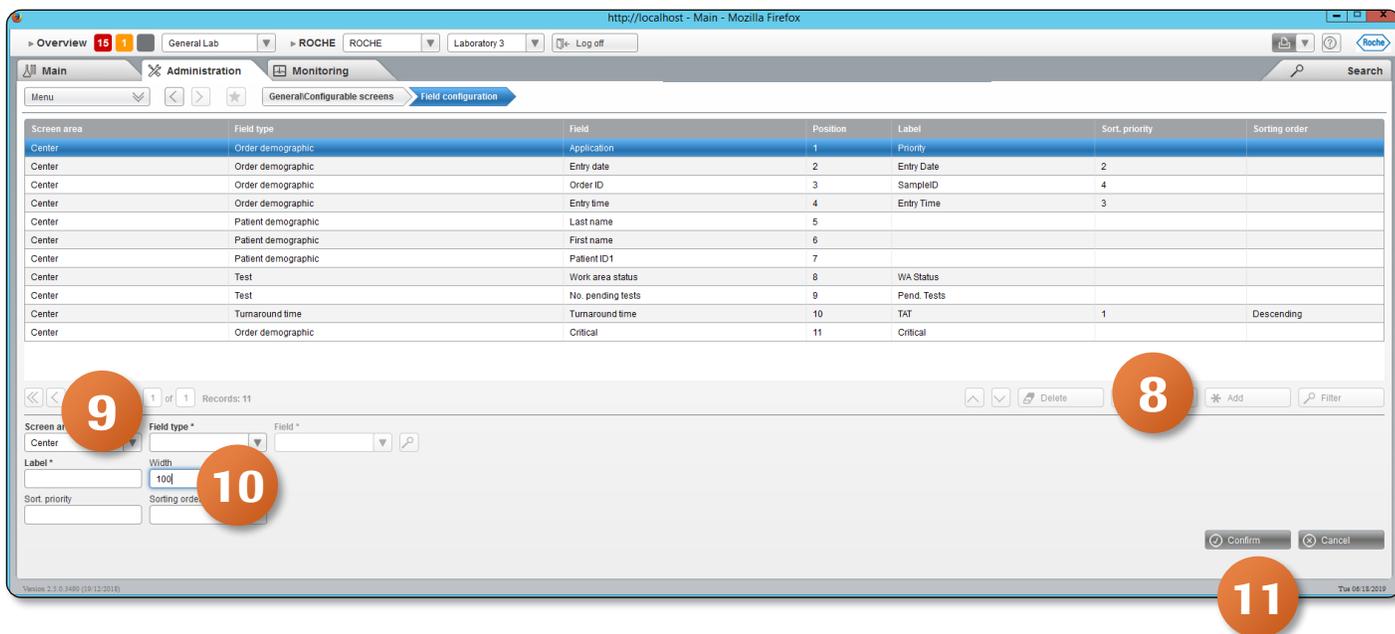

- You also have the option to reorder the location of the columns. To do this, select a row
- Use the **Up and Down** arrows to move to desired location

