



# Searching for a Patient

You can search for patients to edit their record, check their details, patient ID, order information, or results.

1. Once logged into the **cobas® infinity** software, select the **Main** tab
2. Click on **Menu**
3. Navigate to **Queries**
4. Select **Patient search**

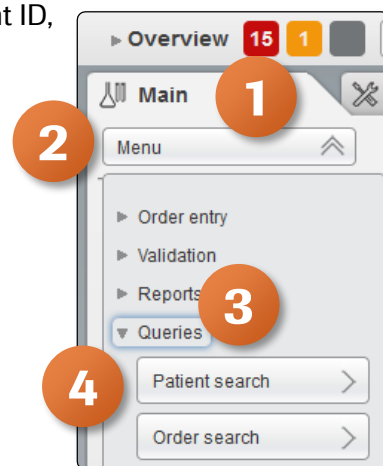

5. Enter the patient demographics you have available, If you don't have any information, you can enter the asterisk (\*) which is known as a wild card. The asterisk (\*) alone will bring up all patients. The asterisk (\*) after a few characters will bring up everything that starts with what was entered. The asterisk (\*) before a few characters will bring up everything that ends with what was entered. The asterisk (\*) before and after the characters will bring up everything that contains what was entered.
6. To search for all patients whose first or last name match the entered patient demographics, select the check box above the **Last name** field
7. Click on **Apply**

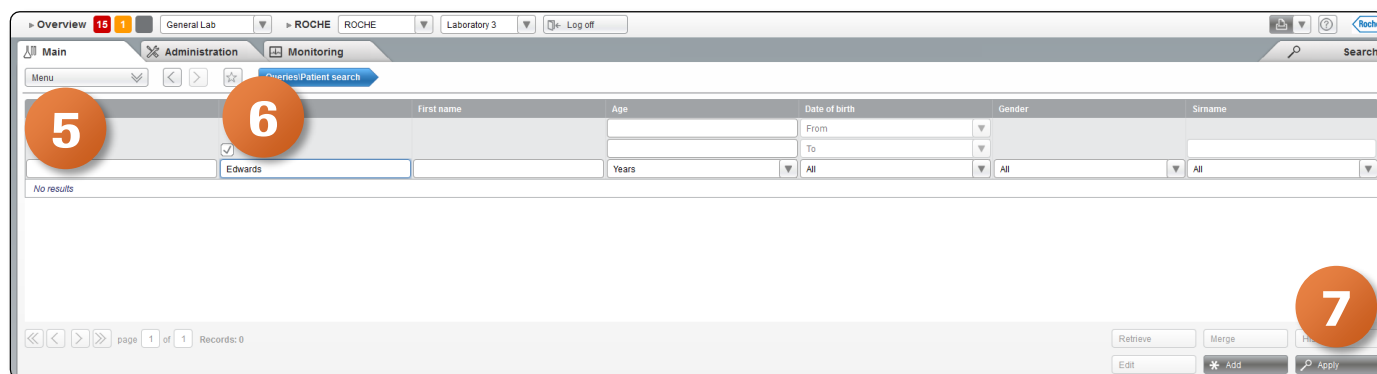

First name	Age	Date of birth	Gender	Surname
Edwards	Years	All	All	All

No results

page 1 of 1 Records: 0

Retrieve Merge Edit Add Apply

# Searching for a Patient

8. Choose the desired patient from top menu
9. The bottom part of the screen displays the orders and pre-orders information belonging to that patient
10. To view an order from that patient, select the order
11. Click on the **Results** button

The screenshot shows the Roche Cobas INFINITY software interface. At the top, there's a navigation bar with 'Overview', 'Administration', and 'Monitoring' tabs. Below this is a search bar and a table of patient records. A red circle with the number '8' highlights the patient selection area. Below the patient table is a table of orders. A red circle with the number '9' highlights the order table. A red circle with the number '10' highlights a specific order row. A red circle with the number '11' highlights the 'Results' button at the bottom right of the order table.

Patient ID1	*Last name	First name	Age	Date of birth	Gender	Surname
123456	Edwards	Theo	57	10/05/1951	Male	

Local dat...	Order ID	Physician	Diagnosis	Service	Destination	Notes	Ward	Location	Critical	Date collected	Time collected	Order status
<input checked="" type="checkbox"/>	Z40000001		Multiple Myeloma					Laboratory 3	Yes			<input checked="" type="checkbox"/>
<input checked="" type="checkbox"/>	Z20000002							Laboratory 3	Yes			<input type="checkbox"/>
<input checked="" type="checkbox"/>	Z30000001							Laboratory 3				<input type="checkbox"/>
<input checked="" type="checkbox"/>	Z20000001							Laboratory 3				<input type="checkbox"/>

12. Depending on your rights, you can access the order in edit or read-only view

The screenshot shows the 'Manual results' view for a specific order. The top section displays 'Order details' and 'Patient details'. Below this is a table of test results. A red circle with the number '12' highlights the 'Results' button at the bottom right of the test results table.

Test	Result	Result alarms	Status	C	Flags	Ref. ranges
BUN	24				0	8 - 25
CA	7.5				0	8.0 - 11.0
CO2	23				0	22 - 29
H	3				0	
I	27				0	
L	11				0	
LACT	4.20				0	0.50 - 2.20
PCT	2.50				0	0.00 - 0.08
GLU	138				0	65.00 - 99.00
CR	1				0	0.70 - 1.30
NA	142				0	135.00 - 145.00
K	5.80				0	3.50 - 5.40
CL	101				0	97.00 - 107.00