



Searching for an Order

You can check information and the status of a previously created order to see which tests were performed and which are pending.

1. Once logged into the **cobas[®] infinity** software, select the **Main** tab
2. Click on **Menu**
3. Navigate to **Queries**
4. Select **Order search**

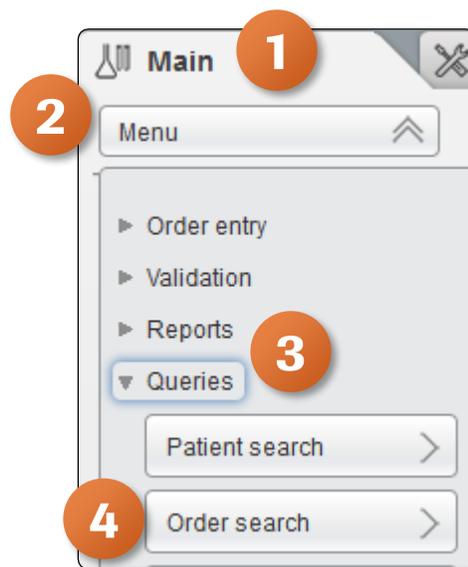


5. From the **Supergroup** drop-down menu, choose the Supergroup the order belongs to
6. Enter the **Year** the order was created. By default the current year shows up. You can edit if necessary
7. Enter either the **Order ID**, **Patient ID**, or **Tube ID**, and then press Enter. If you are unsure of the data, enter an asterisk (*) in any of the fields to search for all information
8. To view the desired order, select it and choose the **Results** button. Depending on your rights, you can access the order in edit or read-only view

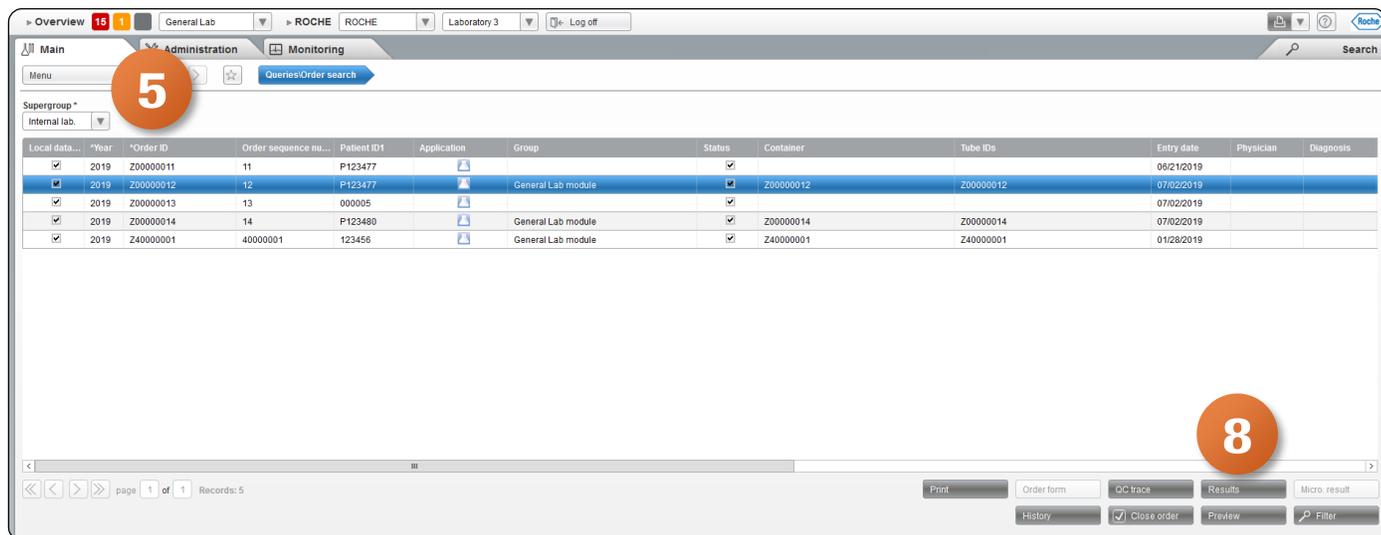

Searching for an Order

9. Navigate back to the **Queries\ Order search** screen by clicking on the bread crumb trail at the top of the screen

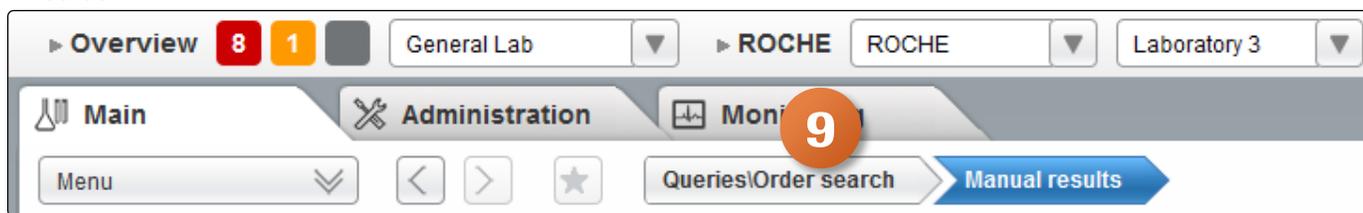

10. Choose the **QC trace** button if you would like to view the QC results related to the tests in the order. This will generate a report with the QC information

