



Setting Up Alarm for Database Size

Setting up the alarm for database size is located in General Parameters. Roche doesn't recommend you change any other General Parameters outside of this alarm.

1. Once logged into the **cobas[®] infinity** software, select the **Administration** tab
2. Click on **Menu**
3. Navigate to **General**
4. Select **General parameters**

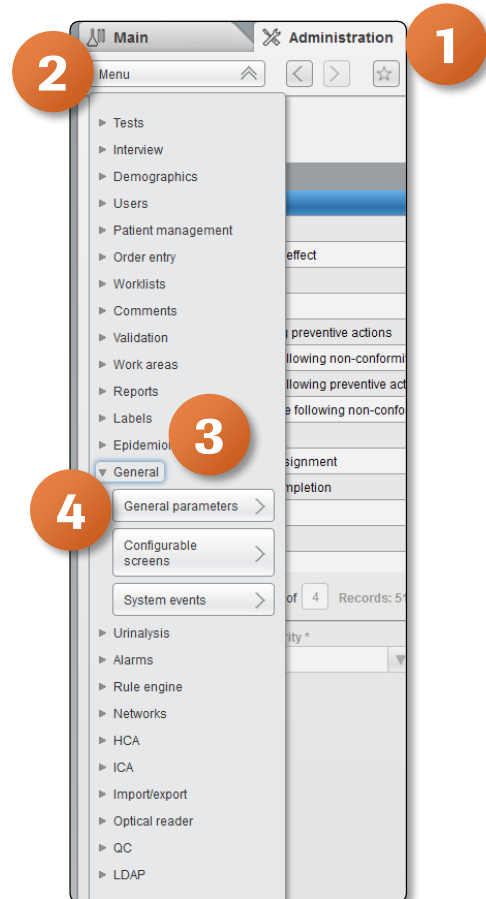


5. Click on **Filter** in the General Parameter screen

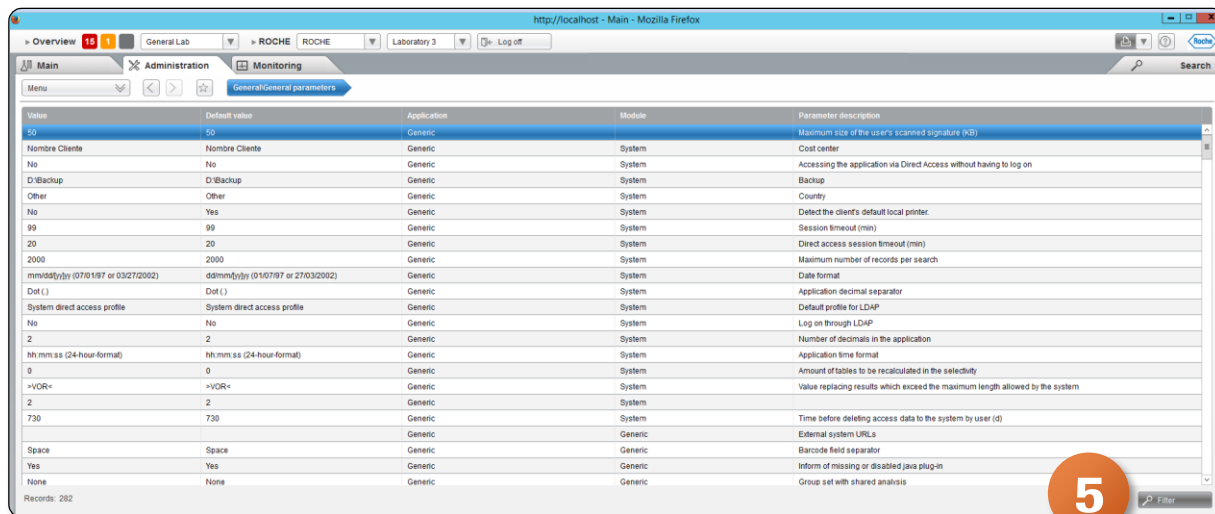

Setting Up Alarm for Database Size

- Place your pointer in the Parameter description box and enter ***alarm***
- Click on **Apply**

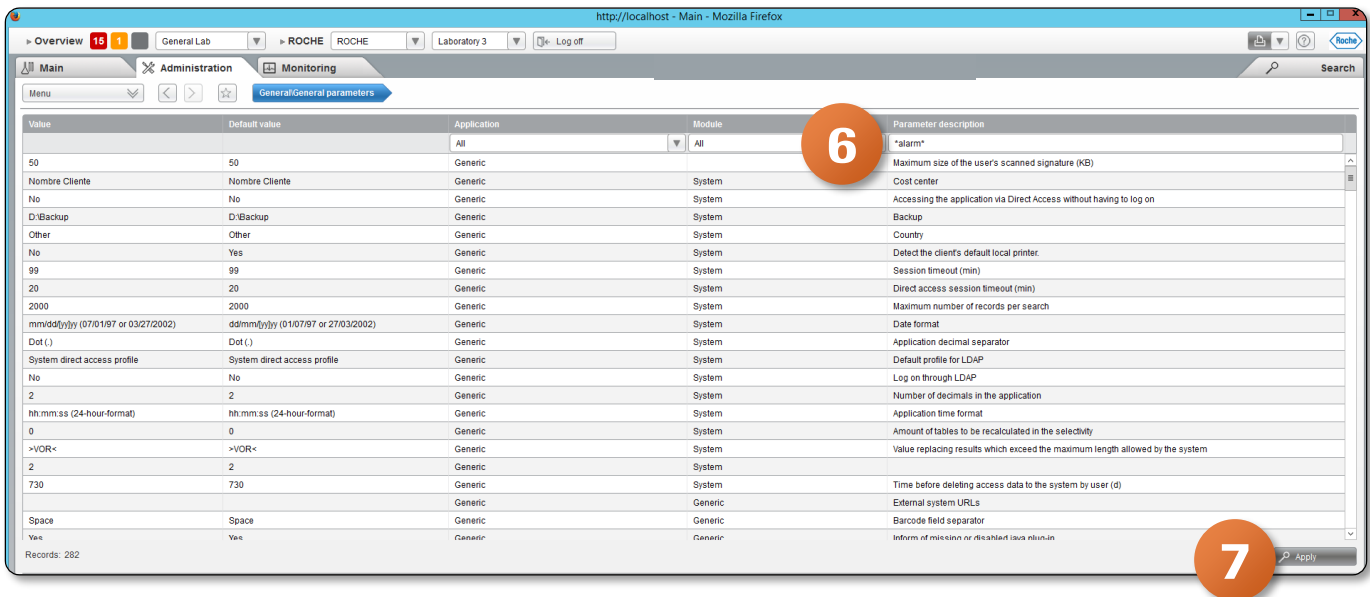

- Double click on the **Percentage of the database size that triggers the maximum size alarm** option in the top portion of the screen
- The bottom portion of the screen now shows the **Value** that will trigger this alarm
- If changes are made, click on **Confirm**

