




Stop and Start Host/LIS Connection(s)

Stop Host/LIS Connection(s)

- Once logged into **cobas[®] infinity** software, select the **Monitoring** tab
- Click on **Menu**
- Navigate to **HCA** (Host Connection Agents)
- Select **Host status**
- A list of configured hosts is shown. The **Enabled** column shows the status of all host connections because of icon
 - The icon indicates the host is waiting to establish the definitive connection. Stop the connection and start again.
- To stop a Host/LIS connection, select the connection from the top of screen
- Click on **Disable**

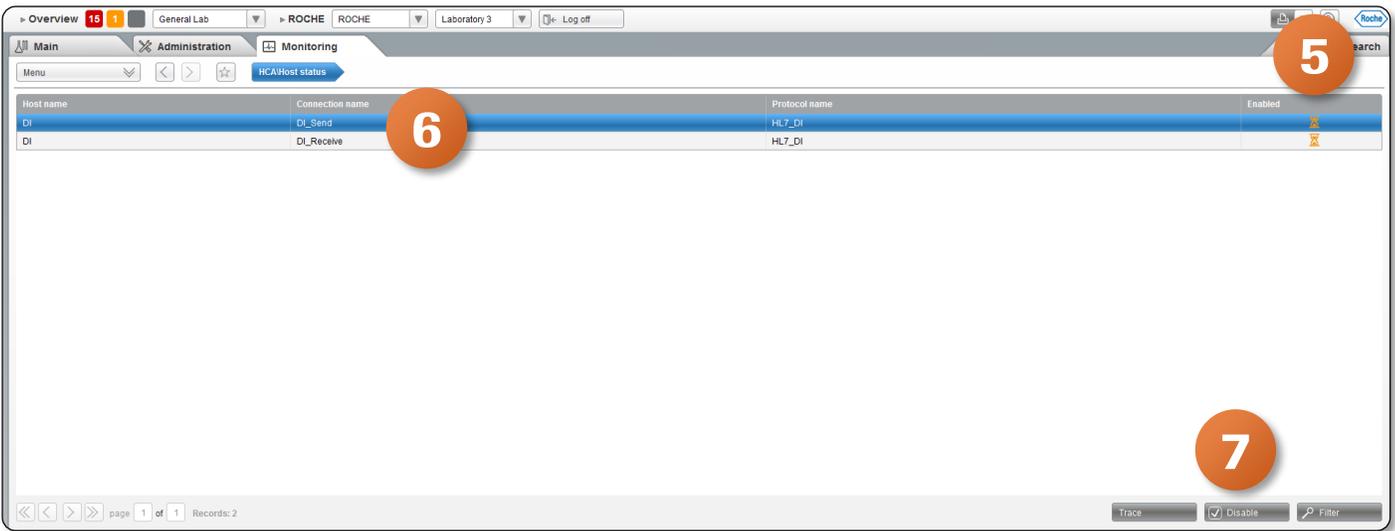


- You will get a pop up asking if you really want to disable
- Click on **Yes**
- You can now see the icon which means this Host/LIS is disabled

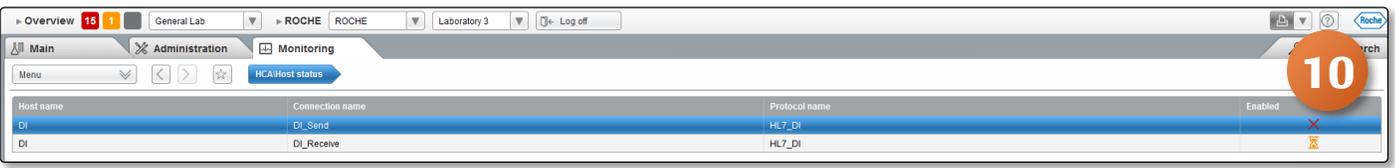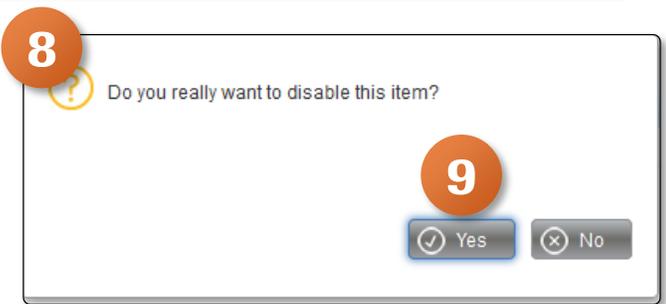

Stop and Start Host/LIS Connection(s)

Start Host/LIS Connection(s)

1. Once logged into **cobas® infinity** software, select the **Monitoring** tab
2. Click on **Menu**
3. Navigate to **HCA** (Host Connection Agents)
4. Select **Host status**
5. A list of configured hosts is shown. The **Enabled** column shows all hosts disabled because of  icon
 - The  icon indicates the host is waiting to establish the definitive connection. Stop the connection and start again.
6. To start a Host/LIS connection, select the connection from top of the screen
7. Click on **Enable**

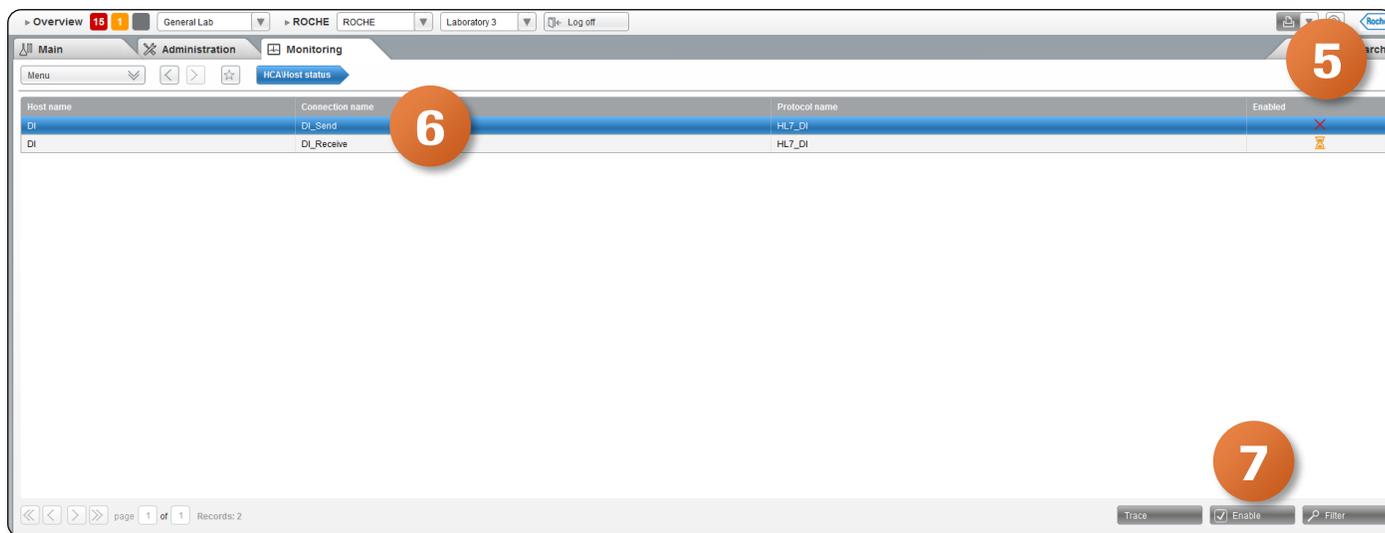

8. You can see that the status is now  showing that the instrument is connected

