
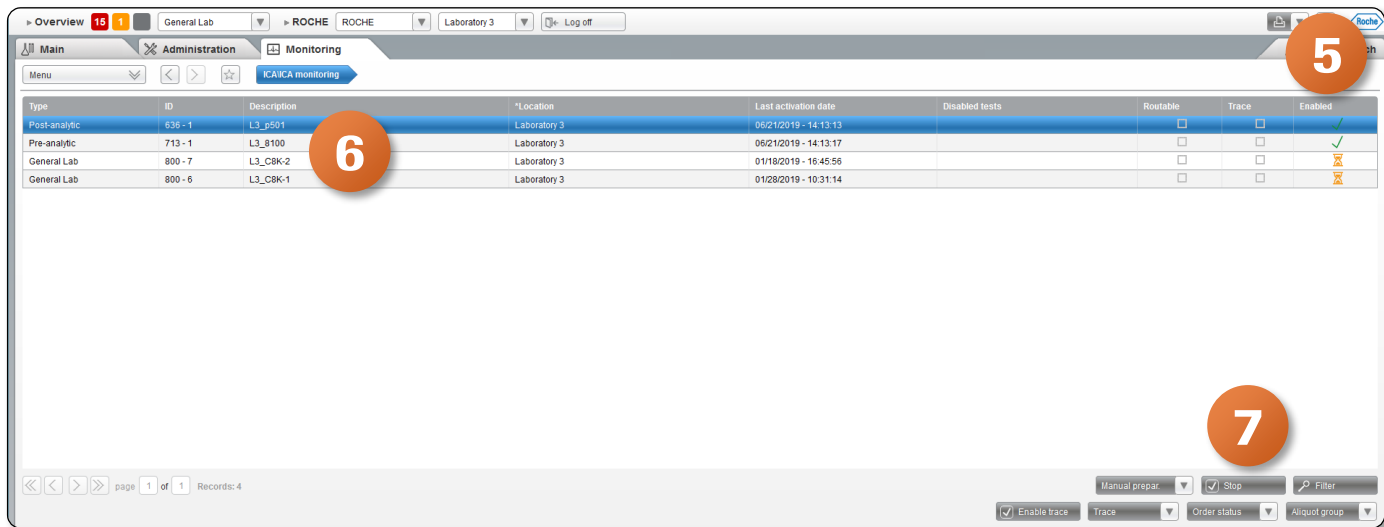




Stop and Start Instrument Connection(s)

Stop Instrument Connection(s)

1. Once logged into **cobas[®] infinity** software, select the **Monitoring** tab
2. Select **Menu**
3. Navigate to **ICA** (Instrument Connection Agents)
4. Select **ICA monitoring**
5. A list of configured instruments is shown. The **Enabled** column shows all instruments connected because of icon
 - The icon indicates the host is waiting to establish a definitive connection. Stop the connection and start again
6. To stop an instrument, select desired instrument from top of screen
7. Click on **Stop**

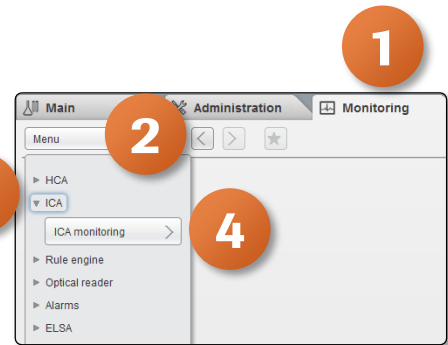


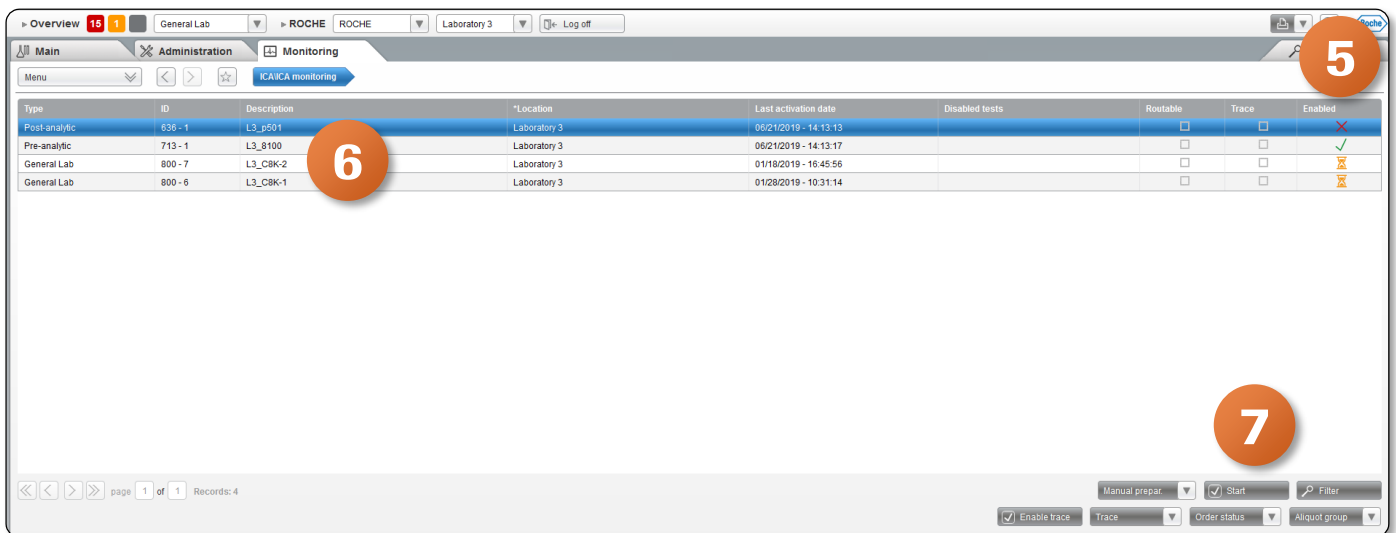
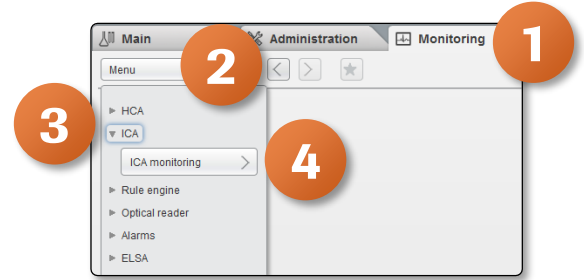
8. You will get a pop up asking if you really want to disable
9. Click on **Yes**
10. You can now see the icon which means this instrument is disabled

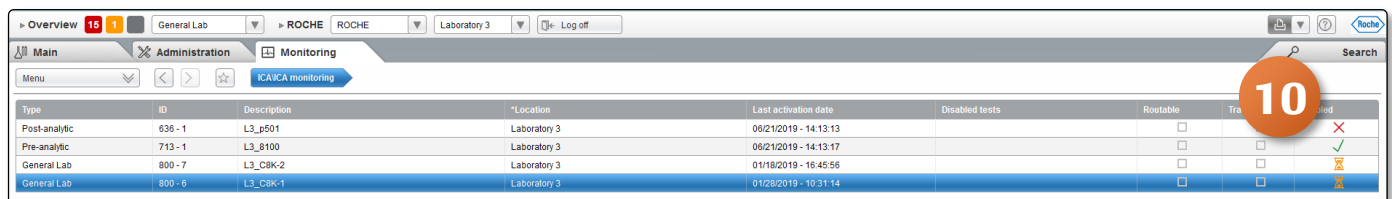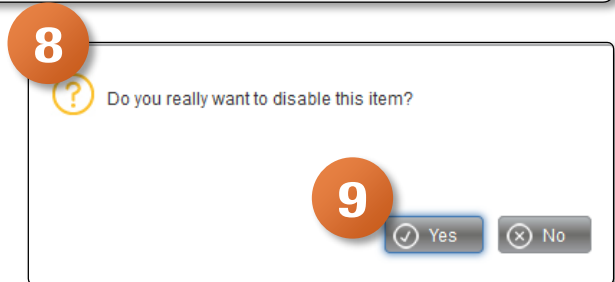

Stop and Start Instrument Connection(s)

Start Instrument Connection(s)

1. Once logged into **cobas® infinity** software, select the **Monitoring** tab
2. Select **Menu**
3. Navigate to **ICA** (Instrument Connection Agents)
4. Select **ICA monitoring**
5. A list of configured instruments is shown. The **Enabled** column shows all disabled instruments because of  icon
6. To start an instrument, select desired instrument from top of screen
7. Click on **Start**



8. You can see that the status is now  showing that the instrument is connected

