

Viewing the Patient's Previous Orders


You can see previous results for a certain patient to check result consistency and give an accurate diagnosis. You can see the patient's previous orders via the Patient history tab

To View the Patient's Previous Orders

1. Once logged into the **cobas[®] infinity** software, select the **Main** tab
2. Click on **Menu**
3. Navigate to **Work areas**
4. Select the Work area you wish to view

Please note: The Work Area names are customized per laboratory. What you see in our example won't necessarily reflect the names of your Work Areas. The above navigation will help you see the names of all Work areas you have created.

5. Double click the desired order from the Monitoring screen to bring up the Validation screen.

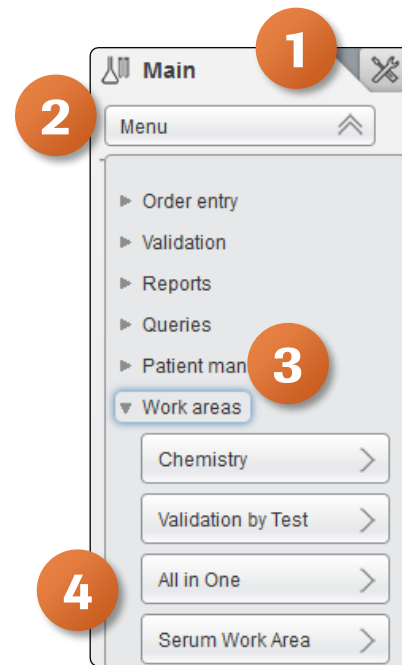

Priority	Entry Date	SampleID	Entry Time	First name	Last name	Patient ID1	WA Status
	01/28/2019	Z40000001	10:38:51	Theo	Edwards	123456	■ ■ ■ ■ ■ ■
	06/21/2019	Z00000011	19:41:42	Brennan	Matthews	P123477	■ ■ ■ ■ ■ ■
	07/02/2019	Z00000012	09:30:41	Brennan	Matthews	P123477	■ ■ ■ ■ ■ ■
	07/02/2019	Z00000014	09:39:01	Estrella	Douglas	P123480	■ ■ ■ ■ ■ ■
	07/11/2019	Z00000015	09:31:54	Vern	Smith	000008	■ ■ ■ ■ ■ ■
	07/11/2019	Z00000016	10:54:45	Adam	Smith	000009	■ ■ ■ ■ ■ ■

Viewing the Patient's Previous Orders

6. Select up to two test check boxes to see their results
7. Choose the **Previous Results** tab
8. Optionally, to see the results from oldest to most recent orders, select the **Show chart from oldest to most recent orders**

The screenshot displays the COBAS INFINITY software interface. On the left, a table lists various test results for a patient. On the right, a chart shows the 'Previous Results' for a specific test (CA). A table below the chart provides details for the two most recent orders.

Test name	Test result	Primary unit	T	Ref. ranges	Critical range	Result alarms	Status	Test alarms
Bicarbonate	7	mmol/L		22 - 29				0
BUN	24	mg/dl		8 - 25				0
Calcium	7.5	mg/dL		8.0 - 11.0				0
Hemolytic	3							0
Icteric	27							0
Lactate	4.20	mmol/l		0.50 - 2.20	> 4.00	▲▲		0
Lipemic	11							0
Procalcitonin	2.50	ng/mL		0.00 - 0.08	> 1.00	▲▲		0
GLUCOSE	138	mg/dL		65.00 - 99.00	< 30.00, > 500.0	▲		0
CREATININE	1	mg/dl		0.70 - 1.30	> 10.00			0
SODIUM	142			135.00 - 145.00	< 125.00, > 155			0
POTASSIUM	5.80			3.50 - 5.40	< 2.50, > 6.50	▲		0
CHLORIDE	101			97.00 - 107.00	< 90.00, > 120.0			0

Test	Result	C	Alarm	Creation date and time	Order ID
CA	9.5			07/24/2018 22:01:16	Z3000002
CA	9.5			07/24/2018 17:25:18	Z3000001